

## **PYROSTAR™ ES-F**

*Multi Test Vial (2.0 mL and 5.2 mL) with Control Standard Endotoxin (CSE)*

**Intended Use:** *Limulus* amebocyte lysate (LAL) is intended for the detection of Gram-negative bacterial endotoxins. PYROSTAR™ ES-F is intended for the qualitative detection of endotoxins by the gel-clot method or quantitative detection by the kinetic turbidimetric method. The quantitative range for the Kinetic Turbidimetric Assay (KTA) is based on the lysate gel-clot labeled sensitivity, refer to the table below:

<b>Gel-clot Labeled Lysate Sensitivity (EU/mL)</b>	<b>KTA Quantitative Range (EU/mL)</b>
0.015	0.001 to 10
0.03 to 0.25	0.01 to 10

### **SUMMARY AND GENERAL INFORMATION**

Endotoxin (lipopolysaccharide or LPS) is a component of the outer membrane of Gram-negative bacteria. Since endotoxins, when injected or implanted, may cause fever and/or shock, the detection of bacterial endotoxins in pharmaceuticals and medical devices is critical.

The LAL assay is the most sensitive method for the detection of bacterial endotoxins currently approved by the U.S. Food and Drug Administration (FDA).<sup>2</sup> The first methodology used for the LAL test was the formation of a gel-clot in the bottom of a glass reaction tube. It has also been observed that the test solution becomes turbid prior to gel-formation. The time required to produce a specified level of turbidity is inversely proportional to the amount of endotoxin in a sample.<sup>1</sup>

A photometric instrument such as the Toxinometer® (FUJIFILM Wako Pure Chemical Corporation) is used to measure the rate of turbidity change. This quantitative measurement procedure is often referred to as the Kinetic Turbidimetric Assay (KTA).

By utilizing these properties, FUJIFILM Wako Chemicals U.S.A. Corporation (FUJIFILM Wako) has developed a LAL assay that can be used either as a quantitative turbidimetric assay or as a qualitative gel-clot assay.

The USP <85> Bacterial Endotoxins Test<sup>6</sup> provides standardized procedures for validation prior to routine use.

The specificity of LAL; however, is not absolute.<sup>7</sup> It has been reported that LAL reacts not only with endotoxin but also with  $\beta$ -1,3-glucan. Although the cascade system activated by  $\beta$ -1,3-glucan has been shown to be different than the one activated by endotoxin,<sup>8</sup> the end result of gel-clot formation is indistinguishable.

The activation of LAL by glucan in a sample can be prevented by adding a large amount of carboxymethylated curdlan (CMC) to LAL. The presence of large amounts of glucan does not interfere with the quantitation of endotoxin. FUJIFILM Wako first made use of these findings by developing an endotoxin-specific buffer (ES-buffer), which contains high concentrations of CMC. When the ES-buffer is used to reconstitute LAL, the LAL reagent becomes endotoxin-specific. PYROSTAR™ ES-F is a new preparation of LAL in which CM-curdlan is co-lyophilized with LAL. Reconstituting PYROSTAR™ ES-F with LAL Reagent Water results in an endotoxin-specific LAL reagent.

## **HISTORY AND BIOLOGICAL PRINCIPLE**

The *in vitro* detection of bacterial endotoxin was pioneered by Levin and Bang<sup>3</sup>. Their findings showed that the blood of the horseshoe crab, *Limulus polyphemus*, clots in the presence of Gram-negative bacteria. They subsequently reported that all the components required for clot formation could be isolated from the circulating amebocytes found in *Limulus* blood.<sup>3,4</sup>

A guideline was released by the FDA in 1987<sup>5</sup> to inform manufacturers of human drugs and biologicals, animal drugs, and medical devices of procedures that FDA considers necessary to validate the use of LAL as an end-product endotoxin test. The FDA guideline was combined with the USP Bacterial Endotoxins Test in 2000 making the USP <85> Bacterial Endotoxins Test the standard method for U.S. manufacturers. The 1987 FDA Guideline was withdrawn in 2011.

## **WARNINGS AND GENERAL PRECAUTIONS**

PYROSTAR™ ES-F is intended for the *in vitro* detection of Gram-negative bacterial endotoxins. Exercise caution when handling LAL because its toxicity is unknown.

This test is not a diagnostic device and is not to be used to determine endotoxin levels in humans for diagnostic purposes.

## **REAGENTS PROVIDED**

**PYROSTAR™ ES-F Reagent:** PYROSTAR™ ES-F is a lyophilized reagent containing *Limulus* amebocyte lysate, buffers, carboxy-methyl curdlan, monovalent and divalent cations.

**Preparation:** Tap the vial on a level surface to ensure that all powder is at the bottom of the vial. Gently remove the stopper and add 2.0 mL or 5.2 mL of LAL Reagent Water (LRW) to the vial. Replace the stopper and swirl gently to dissolve the contents without contacting the stopper with the LAL solution.

**Storage:** Store lyophilized product at 2-10°C. Refrigerate reconstituted reagent at 2-8°C, maximum 8 hours or at -15 ± 5°C for a maximum of 14 days. Reconstituted reagent can only be frozen and thawed once.

**Note:** For sensitivity 0.015 EU/mL reagent, guaranteed stability is for a maximum 4 hours after reconstitution.

**Control Standard Endotoxin (CSE):** A lyophilized reagent that contains refined endotoxin from *E. coli* and is used to confirm LAL reagent sensitivity, validate product test methods, and prepare inhibition or enhancement testing. The CSE contains 500 ng of endotoxin and is tested in comparison to the U.S. Reference Standard Endotoxin (RSE) to determine the conversion factor (EU/ng)

**Preparation:** Release the vacuum by gently pulling up and removing the rubber stopper from the vial. Aseptically add the appropriate volume of LRW to yield a 1,000 EU/mL solution (information located on the Certificate of Analysis). Replace the rubber stopper, invert the vial several times and vortex vigorously for 2 minutes.

**Storage:** Reconstituted CSE can be stored at 2-10°C for 1 month. CSE should not be stored at temperatures below freezing. When the stored solution is used, vigorously vortex the solution for 1 minute prior to use.

#### **MATERIALS AND EQUIPMENT NOT PROVIDED**

**Reference Standard Endotoxin (RSE):** A U.S. Pharmacopeia reference standard defined potency of 10,000 USP Endotoxin Units (EU).

**LAL Reagent Water (LRW):** Endotoxin-free water.

**Endotoxin free consumables:**

- Pipettes tips
- Dilution tubes

**Toxinometer® Method**

- Toxinometer® (FUJIFILM Wako Pure Chemical Corporation)

- Endotoxin-free test tubes and aluminum caps for Toxinometer®

**Gel-clot Techniques**

- Water bath or heating block incubator capable of maintaining  $37 \pm 1^\circ\text{C}$

- Endotoxin-free test tubes and aluminum caps

#### **PREPARATION OF CONTROL STANDARD ENDOTOXIN DILUTIONS**

**Control Standard Endotoxin Preparation**

Based on the information provided on the Certificate of Analysis, prepare a CSE standard to a concentration of 1,000 EU/mL, as described above.

**Note:** If the Pyrostar ES-F kit used has an 8-digit product code, the CSE potency (EU/vial) is present on the CSE vial label.

Vortex the 1,000 EU/mL CSE solution for at least 2 minutes at room temperature. Using the 1,000 EU/mL CSE solution, prepare the endotoxin dilution series as shown in Table 1 or 2. Vortex each tube for 30 seconds between dilutions. Dilutions may be prepared in different volumes if the same ratio is maintained.

**Table 1**  
**Sample Endotoxin Dilution Scheme**  
**(Two-fold dilution series)**

Initial endotoxin (EU/mL)	Volume of LRW added (mL)	Volume of endotoxin added (mL)	Final endotoxin (EU/mL)
1000 (CSE)	3.6	0.4	100
100	3.6	0.4	10
10	3.6	0.4	1
1	2.0	2.0	0.5
0.5	2.0	2.0	0.25
0.25	2.0	2.0	0.125
0.125	2.0	2.0	0.06
0.06	2.0	2.0	0.03
0.03	2.0	2.0	0.015
0.015	2.0	2.0	0.0078
0.0078	2.0	2.0	0.0039

**Table 2**  
**Sample Endotoxin Dilution Scheme**  
**(Ten-fold dilution series)**

Initial endotoxin (EU/mL)	Volume of LRW added (mL)	Volume of endotoxin added (mL)	Final endotoxin (EU/mL)
1000 (CSE)	3.6	0.4	100
100	3.6	0.4	10
10	3.6	0.4	1
1	3.6	0.4	0.1
0.1	3.6	0.4	0.01
0.01	3.6	0.4	0.001

### **SPECIMEN COLLECTION AND PREPARATION**

LAL reactions are pH sensitive, requiring that the LAL and sample mixture have a pH of 6.0 to 8.0. PYROSTAR™ ES-F contains buffering components that help bring the test mixture within the pH range in most cases. If a pH adjustment is necessary, the pH of the sample should be adjusted with endotoxin-free HCl or NaOH.

### **PRODUCT INTERFERENCE**

Prior to using the LAL test for the routine release of products, it is necessary to validate the absence of product interference by performing an inhibition/enhancement test for each product type. A product is determined to be non-interfering by assaying a sample of the product spiked with a known amount of endotoxin and detecting 50 – 200% recovery of the spiked endotoxin.

**Kinetic Turbidimetric Assay (KTA):** Prepare a standard curve that covers the testing range. Spike the product to a concentration of endotoxin that is equal to or near the middle of the standard curve. The product is determined to be non-interfering if the percent recovery reported is 50 to 200% of the spiked endotoxin concentration. For example, if the standard curve is 0.1 EU/mL, 1.0 EU/mL 10 EU/mL, the product should be spiked with endotoxin to yield a concentration of 1.0 EU/mL. The acceptable endotoxin recovery is between 0.5 and 2.0 EU/mL.

**Gel-clot Assay:** Two independent CSE dilution series are prepared using the product and LRW as diluents, respectively. Assay the CSE dilution series prepared in LRW in duplicate (n=2), and the CSE dilution series prepared in product in quadruplicate (n=4). The geometric mean endpoints of both dilution series must fall within two-fold of the labeled endotoxin sensitivity.

## **KINETIC TURBIDIMETRIC ASSAY PROCEDURE**

A Kinetic Turbidimetric Assay (KTA) can be performed using the Toxinometer® and Toximaster software.

In addition to the product being tested, a valid assay will include endotoxin standards bracketing the analysis range, positive product controls, and negative controls. All test values need to be determined from at least duplicate samples.

**Toxinometer® Method:** Aseptically transfer 0.1 mL of PYROSTAR™ ES-F reagent into each endotoxin-free Toxinometer® tube. Add 0.1 mL of product sample or control to the respective test tube and quickly vortex (~3 seconds) the contents without foaming. Place the tube into the Toxinometer® and proceed to the next Toxinometer® tube. After all reaction tubes are added to the Toxinometer®, the Toximaster software will calculate the endotoxin concentration for each sample and generate the regression analysis based on the data for standard curve.

See "<85> Bacterial Endotoxins Test" in The U.S. Pharmacopeia and refer to sections describing standard curve and preparation of solutions.<sup>6</sup>

### **Calculation of endotoxin concentration**

During a KTA reaction the turbidity of the test solutions is continually monitored by the Toxinometer®. The time required for a sample to reach a predetermined threshold absorbance level over background is measured. This time is referred to by the Toximaster® software as Ta (activation time).

The software produces a x versus y axis correlation of the Ta of each standard (y axis) with its corresponding endotoxin concentration (x axis). An example of a standard series and recovery of an endotoxin spike of 1.0 EU/mL in product is presented below:

#### ***Representative Analysis***

<b>Standards</b>	<b>Endotoxin concentration (EU/mL)</b>	<b>Ta (min)</b>	<b>Log endotoxin Concentration</b>	<b>Log Ta (min)</b>
NC	0	NR	-	-
STD1	10.0	8.0	1.000	0.903
STD2	1.0	13.2	0.000	1.121
STD3	0.1	23.6	-1.000	1.373

slope	-0.2349
y-intercept	1.132
coefficient of correlation	-0.999

		Ta (min)	Log Ta (min)	Calculated EU/mL	% Recovery
Product 1	NPC1	NR	-	-	-
	PPC1	13.6	1.134	0.981	98.1%
Product 2	NPC2	NR	-	-	-
	PPC2	13.0	1.114	1.193	119.3%

NR= Nonreactive

In this example, the positive product control (PPC) for each product sample yielded an endotoxin recovery that was consistent with the spiked level indicating there were no detectable product enhancement or inhibition. The negative product control (NPC) and negative control (NC) showed significantly lower endotoxin levels than the lowest standard endotoxin concentration.

### Performance characteristics

The linearity of the standard curve used to determine endotoxin with the concentration range must be verified. At least three endotoxin standards spanning the desired concentration range and a LRW water blank should be assayed in triplicate. Refer to the Assurance of Criteria for the Standard Curve in USP. The absolute value of the coefficient of correlation,  $r$ , must be greater than or equal to 0.980.<sup>6</sup>

### GEL-CLOT ASSAY PROCEDURE:

To ensure both the precision and accuracy of the PYROSTAR™ ES-F test, the sensitivity indicated on the label must be verified as described in the USP <85> Bacterial Endotoxins Test.<sup>6</sup>

### Test for Confirmation of Labeled Sensitivity

The sensitivity indicated on the PYROSTAR™ ES-F label can be confirmed by preparing a two-fold dilution series of RSE or CSE that brackets the stated sensitivity. The dilution series needs to be assayed in quadruplicate and include negative controls. The sensitivity is then determined by calculating the geometric mean of the obtained endpoints. An example of PYROSTAR™ ES-F with a labeled sensitivity ( $\lambda$ ) of 0.125 EU/mL is described below:

PYROSTAR™ ES-F ( $\lambda$ ) = 0.125 EU/mL

Replicate	0.25	0.125	0.06	0.03	NC	Endpoint	Log <sub>10</sub> Endpoint
1	+	+	+	-	-	0.06	-1.222
2	+	+	-	-	-	0.125	-0.903
3	+	+	-	-	-	0.125	-0.903
4	+	+	+	-	-	0.06	-1.222
Mean (Log <sub>10</sub> Endpoint)							<b>-1.062</b>
Antilog <sub>10</sub> (Mean Log <sub>10</sub> Endpoint)							<b>0.087</b>

Convert each endpoint of the quadruplicate assay to its  $\log_{10}$  value. The individual  $\log_{10}$  endpoint values are averaged. The geometric mean is  $\text{antilog}_{10}$  of the mean. The mean  $\log_{10}$  endpoint value in the above example is -1.062. The geometric mean is 0.087 EU/mL, which confirms the labeled sensitivity is between  $0.5\lambda$ - $2\lambda$  (0.06-0.25 EU/mL) and all negative controls are negative.

### Routine Testing with PYROSTAR™ ES-F

Each assay should include a negative control (NC),  $2\lambda$  positive control ( $2\lambda$  PC),  $2\lambda$  positive product control ( $2\lambda$  PPC) and the product to be tested all in duplicate, where  $\lambda$  is the PYROSTAR™ ES-F sensitivity.

1. Add 0.1 mL of the reconstituted PYROSTAR™ ES-F to each assay tube.
2. Aseptically transfer 0.1 mL of each product sample or control into each assay tube, beginning with the negative control and ending with the highest endotoxin concentration.
3. Quickly mix the contents of the tubes without foaming and incubate them undisturbed in a  $37 \pm 1^\circ\text{C}$  heating block or heated water bath for 60 minutes  $\pm 2$  minutes.
4. After incubation, examine each tube for gelation. Each tube in the gel-clot method is interpreted as either positive or negative. A positive test is defined as the formation of a firm gel capable of maintaining its integrity when the assay tube is inverted  $180^\circ$ . A negative test is characterized by the absence of gel or by the formation of a viscous mass which does not hold when the assay tube is inverted.
5. Test results are only valid when the  $2\lambda$  PC and  $2\lambda$  PPC are positive, and the NC is negative.

### Determination of endotoxin in an unknown sample

To determine the endotoxin concentration of an unknown solution, test two-fold dilutions of the product until an endpoint has been reached. Calculate the geometric mean dilution factor and multiply by the labeled endotoxin sensitivity.

PYROSTAR™ ES-F ( $\lambda$ ) = 0.125 EU/mL

Product Dilution	1/2	1/4	1/8	1/16	1/32	1/64	
Dilution Factor	2	4	8	16	32	64	$\text{Log}_{10}(\text{endpoint dilution factor})$
Replicate 1	+	+	-	-	-	-	0.602
Replicate 2	+	+	+	-	-	-	0.903

Endpoint Dilution Factor (DF)	$\text{Log}_{10}$ Endpoint DF
4	0.602
8	0.903
Mean ( $\text{Log}_{10}$ Endpoint DF)	0.753
$\text{Antilog}_{10}$ (Mean $\text{Log}_{10}$ endpoint DF)	5.66

Endotoxin Concentration (EU/mL) =  $0.125 \text{ EU/mL} \times 5.66 = 0.707 \text{ EU/mL}$  ✓

## **BIBLIOGRAPHY**

1. Cooper, J.F., Levin, J., and Wagner, H.N. "Quantitative Comparison of in vitro and in vivo Methods for Detection of Endotoxin" *Journal of Laboratory and Clinical Medicine*, 78, p.128 (1971).
2. Hochstein, H.D. "The LAL Test versus the Rabbit Pyrogen Test for Endotoxin Detection: Update '87." *Pharm. Technol.*, 11(6), p. 124 (1987).
3. Levin, J. and Bang, F.B. "Clottable protein in Limulus: Its Localization and Kinetics of its Coagulation by Endotoxin." *Thrombosis et diathesis haemorrhagica.*, 19, p. 186 (1966).
4. Tai, J.Y. and Liu, T.Y. "Studies on Limulus Lysate, Isolation of Pro-clotting Enzyme." *Journal of Biological Chemistry*, 252, p.2176 (1977).
5. Guideline on Validation of the Limulus Amebocyte Lysate Test as an End-Product Endotoxin Test of Human and Animal Parenteral Drugs, Biological Products and Medical Devices. U.S. Food and Drug Administration, 1987.
6. "<85> Bacterial Endotoxins Test." The United States Pharmacopeia, most current version. U.S. Pharmacopeial Convention, Rockville, MD.
7. Cooper, J.F., Weary, M.E., and Jordon, J.T. "The Impact of Non-endotoxin LAL-Reactive Materials on *Limulus* Amebocyte Lysate Analysis." *PDA J Pharm Sci Technol.*, 51(1), (1997).
8. Tsuchiya, M., Oishi, H., Takaoka, A., Fusamoto, M. and Matsuura, S. "Discrimination between endotoxin and (1 – 3) beta-D-glucan using turbidimetric kinetic assay with *Limulus* amebocyte lysate." *Chem Pharm Bull* (Tokyo), 38(9). p. 2523 (1990).

### **Manufactured by:**

FUJIFILM Wako Chemicals U.S.A. Corporation  
LAL Division  
1600 Bellwood Road  
Richmond, VA 23237

US License No. 2093

### **For Customer Complaints, Contact:**

FUJIFILM Wako Chemicals U.S.A. Corporation  
Quality Assurance in LAL Division  
+1-804-271-7677